

Art. No. 88-52000

- DE** Bedienungsanleitung
- GB** Operating Instructions
- FR** Mode d'emploi
- NL** Handleiding
- IT** Istruzioni per l'uso
- ES** Instrucciones de uso
- PT** Manual de utilização

| | | |
|-----------|------------------------------------|-----------|
| DE | Bedienungsanleitung..... | 4 |
| GB | Operating Instructions..... | 8 |
| FR | Mode d'emploi..... | 12 |
| NL | Handleiding..... | 16 |
| IT | Istruzioni per l'uso..... | 20 |
| ES | Instrucciones de uso | 24 |
| PT | Manual de utilização..... | 28 |

Fig. 1

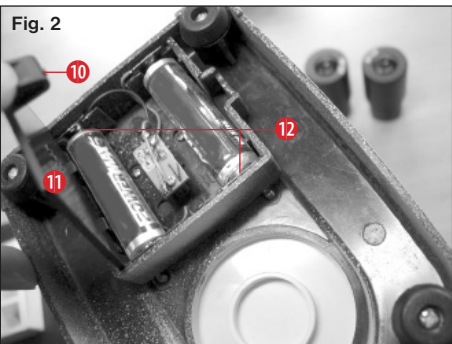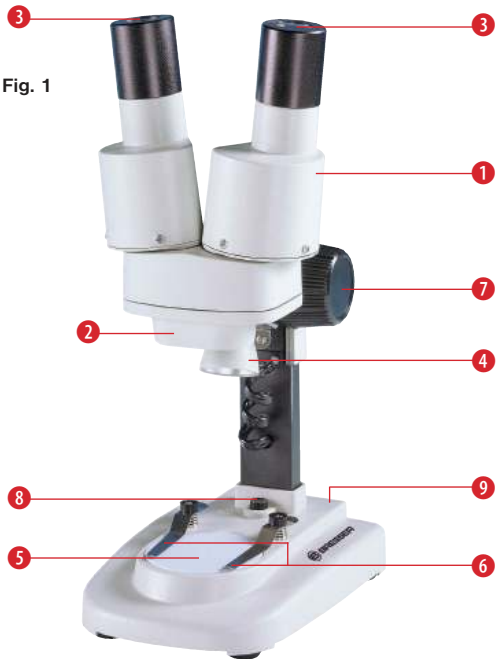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DANGER to your child!

Aids with sharp edges and tips are sometimes used with this device. Please store the device and all of its accessories and aids out of the reach of children. There is a risk of **INJURY**.

This device contains electronic components that are powered by either a mains connection or batteries. Never leave a child unsupervised with this device. The device should only be used as per these instructions otherwise there is a serious **RISK of ELECTRICAL SHOCK**.

Batteries should be kept out of children's reach. When inserting batteries please ensure the polarity is correct. Leaking or damaged batteries can cause injury if they come into contact with the skin. If you need to handle such batteries please wear suitable safety gloves.

Children should only use this device under supervision. Keep packaging materials (plastic bags, rubber bands, etc.) away from children. There is a risk of **SUFFOCATION**.

FIRE-/ DANGER OF EXPLOSION!

Do not expose the device to high temperatures. Use only the mains adapter supplied or those battery types recommended. Never short circuit the device or batteries or throw into a fire. Exposure to high temperatures or misuse of the device can lead to short circuits, fire or even explosion!

RISK of material damage

Never take the device apart. Please consult your dealer if there are any defects. The dealer will contact our service centre and send the device in for repair if needed.

Do not subject the device to temperatures exceeding 60° C.

DISPOSAL

Dispose of the packaging materials properly, according to their type (paper, cardboard, etc). Contact your local waste disposal service or environmental authority for information on the proper disposal.

Do not dispose of electronic devices in the household garbage!

As per the Directive 2002/96/EC of the European Parliament on waste electrical and electronic equipment and its adaptation into German law, used electronic devices must be collected separately and recycled in an environmentally friendly manner.

Empty old batteries must be disposed of at battery collection points by the consumer. You can find out more information about the disposal of devices or batteries produced after 01.06.2006 from your local waste disposal service or environmental authority.

In accordance with the regulations concerning batteries and rechargeable batteries, disposing of them in the normal household waste is explicitly forbidden. Please pay attention to dispose of your used batteries as required by law - at a local collection point or in the retail market (a disposal in domestic waste violates the Battery Directive).

Batteries that contain toxins are marked with a sign and a chemical symbol. „Cd“ = cadmium, „Hg“ = mercury, „Pb“ = lead.

Cd¹

Hg²

Pb³

¹ battery contains cadmium

² battery contains mercury

³ battery contains lead

EEC conformity explanation

CE Meade Instruments Europe GmbH & Co KG, resident in 46414 Rhede/Westf., Gutenbergstr. 2, Germany, explains the agreement with in the following specified EEC guidelines for this product:

EN 61010-1:2001

Product description:

StereoMicro

Model /Description:

Microscope STX

Rhede, 05-05-2009

Meade Instruments Europe GmbH & Co. KG

Helmut Ebbert

Managing director

Your microscope is made up of the following parts (Fig. 1-4):

- 1 Binocular view piece (with eyepiece supports)
- 2 Stereo objective
- 3 Eyepieces
- 4 Illumination (LED)
- 5 Microscope stage (with object plate)
- 6 Stage clips
- 7 Focus knob
- 8 On/off switch
- 9 Battery compartment
- 10 Battery compartment clamp closure
- 11 Battery compartment cover
- 12 Battery contacts (flat sheets and spring terminals)
- 13 Diffusion cap

Location

Before you start, choose an ideal location for using your microscope. It is important that you have enough light available (window, desk lamp).

Simple Observation

Place the object to be observed (for example, a rock) in the middle of the microscope stage (5). When the object is large enough, it is best to fix it in place with the stage clips (6).

Now, take a look through the binocular eye lens (1) and change the distance of the eyepiece (3), so that you have a circular, clear image.

The sharpness is adjusted by slowly turning the focus knob (7).

Illumination

Remove both eyepieces (3) from the eyepiece supports (1) and turn the entire microscope over. In the base of the microscope there is a battery compartment (9). Push the clamp closure (10) and open the battery

compartment by pulling the cover (11) open diagonally.

Place the batteries in the compartment so that the flat minus poles (-) press against the spring terminal and the plus poles (+) are touching the flat contact sheets (12).

Close the battery compartment with the cover (11) and turn the microscope around again. Place the eyepieces back on. You can not turn on the light (4) with the on/off switch (8) on the base of the microscope.

TIPS on cleaning

Remove the device from its energy source before cleaning (remove plug from socket / remove batteries).

Clean the exterior of device with a dry cloth. Do not use cleaning fluids so as to avoid causing damage to electronic components.

Protect the device from dust and moisture. Store the device in the bag supplied or in its original packaging. Batteries should be removed from the device if it is not going to be used for a long period of time.

Clean the lens (objective and eyepiece) only with the cloth supplied or some other soft lint-free cloth (e.g. micro-fibre). Do not use excessive pressure - this may scratch the lens.

Dampen the cleaning cloth with a spectacle cleaning fluid and use it on very dirty lenses.

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